

### **GeneReleaser® REVERSE TRANSCRIPTASE Protocol**

1. Accomplish lysis in a standard PCR tube by one of the two GeneReleaser® protocols (Thermocycle or Microwave Lysis).
2. Centrifuge the tube at max rpm in a microfuge.
3. Take an aliquot of the supernatant from 2 above and perform the Reverse Transcriptase (RT) reaction according to the procedure of your RT vendor.
4. Following RT reaction inactivate RT and take suitable aliquot of the RT reaction and perform PCR according to your standard protocol.

If you have further questions, please contact us at [support@bioventures.com](mailto:support@bioventures.com). You may also contact us via fax (1-877-286-0330) or through our toll-free phone number (1-877-852-7846).