

GeneReleaser® DNA AMPLIFICATION FROM SALIVA Protocol

1. Dilute saliva 1:10 with normal saline
2. Microfuge for 1 minute at ~ 14,000 X g.
3. Discard supernatant.
4. Treat pellet per microwave protocol.
5. Amplify per standard protocols.
6. Serial dilutions of the saliva suspension may be required to establish optimum preparation.

If you have further questions, please contact us at support@bioventures.com. You may also contact us via fax (1-877-286-0330) or through our toll-free phone number (1-877-852-7846).